



This service procedure details the replacement of the LCD Screen Assembly (X90). For further assistance, please contact Technical Services: <https://www.ankom.com/contact/technical-services>


1. Power down the instrument and unplug to ensure safety.
2. Remove the back panel by removing the ten nuts that hold it in place. Set the panel and nuts aside for reassembly.

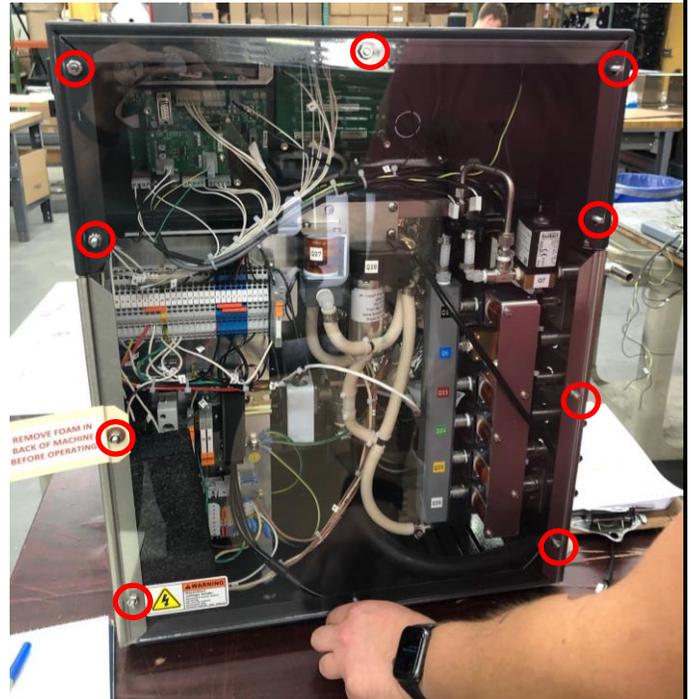


3. Using a 11/32 socket wrench, remove the two screws holding the LCD Screen Assembly in place. Set aside the screws for reassembly.

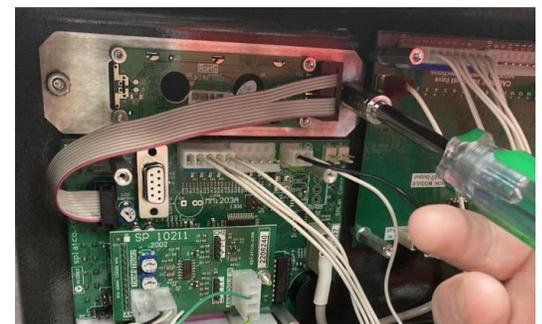

4. Detach the LCD Screen Assembly from the chip board by unclipping the lefthand portion of the connecting wire. Remove the assembly from its position by pulling away gently.

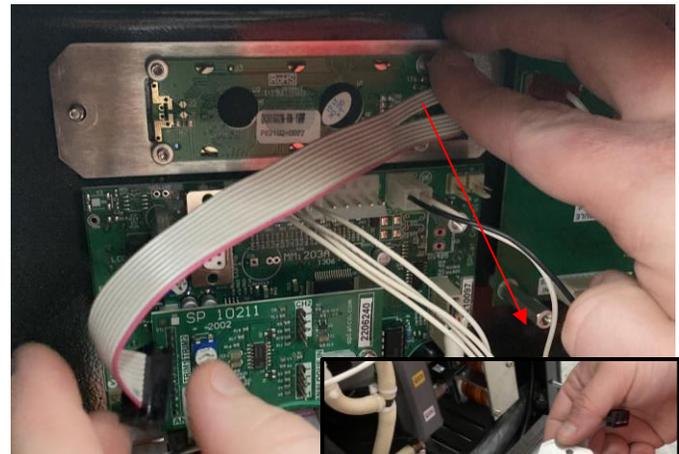


5. Obtain the new LCD Screen Assembly (SKU: #X90).

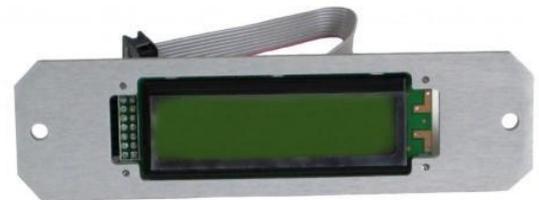

6. Attach the new assembly by holding the top portion in place and reattaching the lefthand portion of the wiring.



7. Using the 11/32 socket wrench, reattach the two screws that hold the assembly in place.

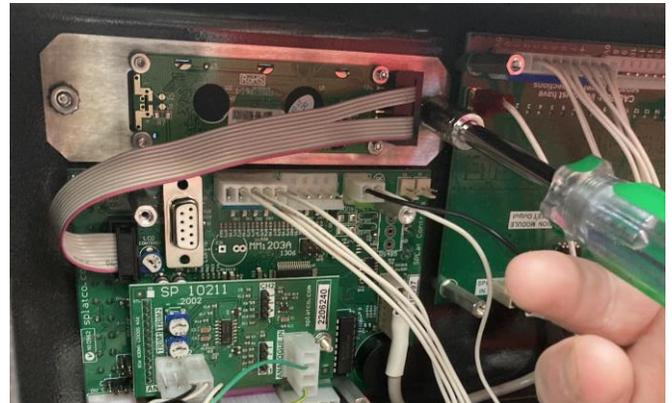

8. Reattach the back panel and install the ten screws back in place.

