Leak Check at Fittings

Service Procedure #003

To check for leaks at the fittings, do the following:

- .) Disconnect the solution supply tubes on the left and right sides and start a run.
- 2) When the screen says "Filling" pour in 2100ml of hot soapy water (dish soap works best) over the bag suspender to cover the level sensor. Agitation will begin. Close the lid.
- 3) Place the open ends of the drain tube and vent line in separate small buckets of water. Do you notice "bubbling" as the instrument heats?
- 4) In the back of the instrument do you notice any bubbles in the A and B Clear Valve Tubing as the instrument heats and all fittings marked A-N?

Check Lid (front of A2000)

(pg 1 of 1) 06/06/2019 e

