


1. Remove the A07 Water Filter Assembly from the 6"x6" plastic bag. This should come with a nut (Red Arrow) already on it.



2. If an A13 Copper Tube Assembly was purchased, the nut (Red Arrow) comes already installed on the Copper tubing. Remove the nut from the A07 Water Filter (This is an extra and can be saved in case the one being used is damaged) and use the one already installed on the Copper tubing Assembly. When connecting it to the A07 Water Filter, Take notice of the direction of the ferrule (Blue Arrow).

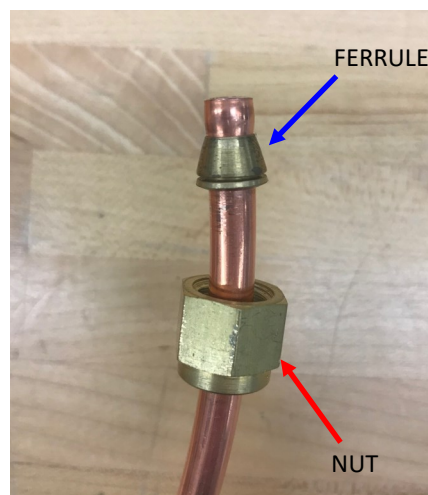

3. Install the Water Filter Assembly (A07) to the port labeled "WATER SUPPLY" on the left side of the DELTA/A2000 Instrument.

