


Controller (Splat) Replacement A2000/A2000I

Caution: Be careful not to discharge static electricity when handling the circuit board. Damage to the program will result. Use a ground strap if possible.

1. Unplug the Instrument.
2. Remove the cabinet back.
3. Spread the small tab on the connector with a small screwdriver or your thumb and disconnect the connectors (1,2,3 and 4,5,6). (Note: Your A2000 may not have a wire in connector # 2. See page 2 for example, installation instructions are the same.)
4. Disconnect the screen and pressure cable (no tab on plug) (7,8).
5. Remove four mounting standoffs.
6. Remove the Splat Controller.
7. Install the new Splat Controller with four standoffs and carefully plug in connectors 1-8. Make sure the connectors are plugged into the same position (match colors).
8. Reattach the cover and cabinet back.

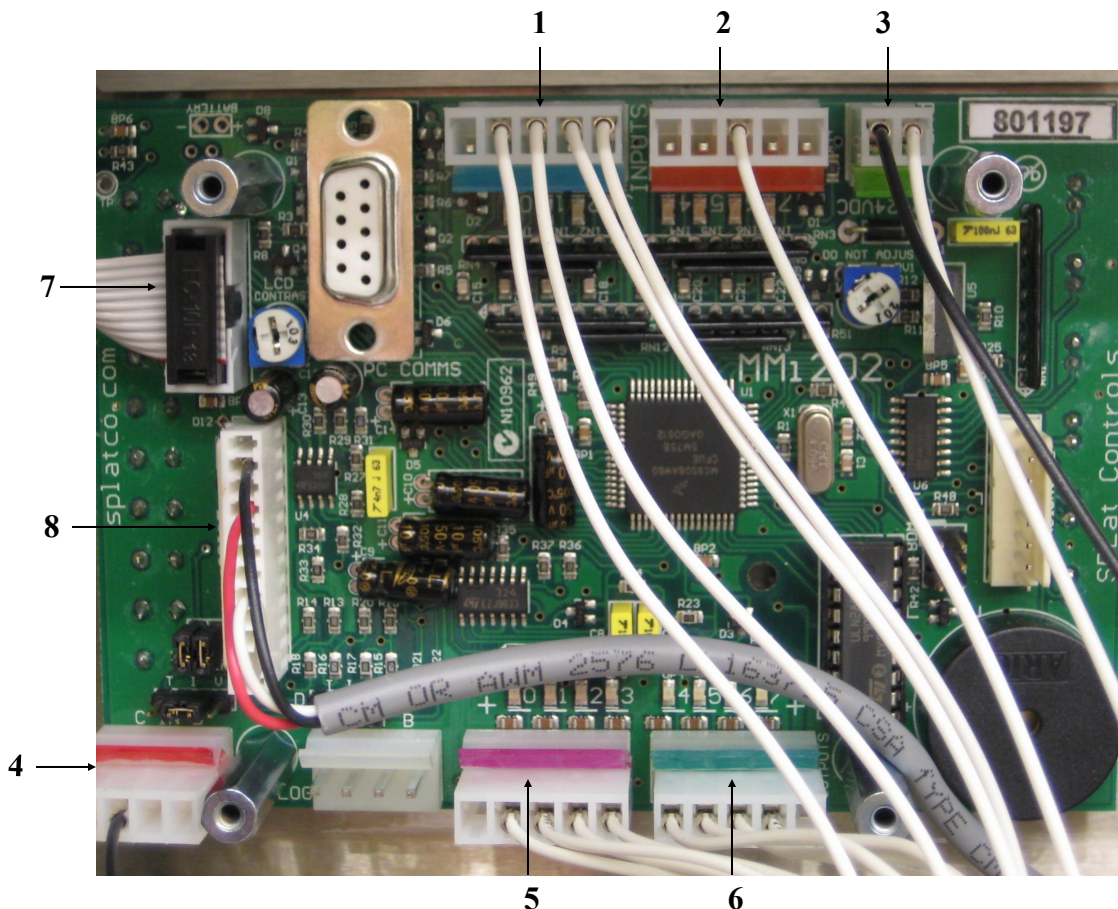

Controller (Splat) Replacement A2000/A2000I

(Note: A2000 instruments produced after 2009 do not have a wire in connector #2. See picture below for example)

