

A01 Cubetainer Shelf Installation A2000/A2000I

Service Procedure **AS024**Rev. C Date: 6/25/19 e

1. Lay the shelf as shown in the picture on right. Remove the six nuts from inside the Shelf Platform (red arrows).



2. Attach the Left and Right Support Brackets, and fasten the nuts back on. Make sure the brackets are oriented correctly.



3. Using a screw driver, remove the screws (red arrows from the Support Post.



4. Secure the Triangular Back Brace to the support post so the outside corners face away from the leveling mount. Replace the screws.





A01 Cubetainer Shelf Installation A2000/A2000I

Service Procedure **AS024**Rev C Date: 6/25/19

5. With the inside facing up, attach the Support Post into the center of the Shelf Platform (red arrows).



6. Remove six nuts from the back of the A2000 (red arrows).



7. Slide on the Left and Right Support Brackets and the Triangular Back Brace as shown (red arrows), then replace the nuts. Adjust the Leveling Foot (blue arrow) as needed to support the bracket. The foam strip should rest on the top of the A2000.







