


Chemical Inlet Port Replacement A2000/A2000I

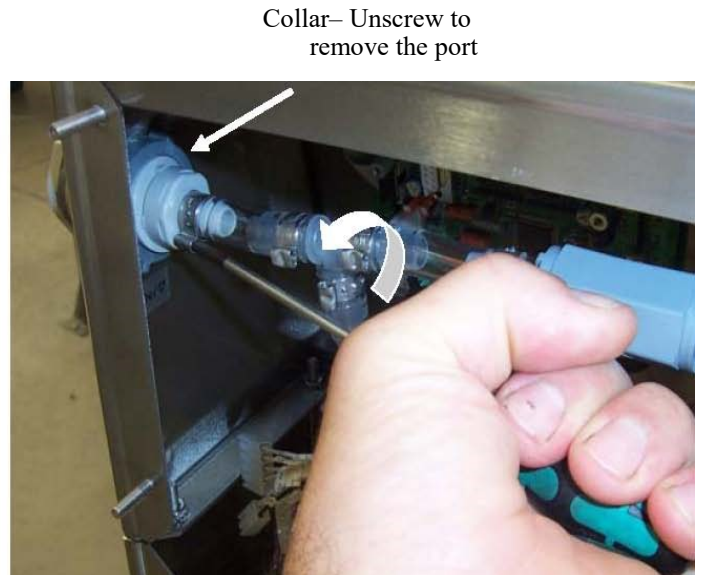

1. Remove the back panel of the A2000.
2. Push a small screwdriver into the crimp of the clamp securing the broken inlet port.
3. Twist the screwdriver as shown to open the crimp.
4. Remove the tube from the port barb and discard the old clamp.
5. Remove the collar holding the inlet port to the cabinet.
6. Install the new inlet port and collar in the cabinet. Slide the screw clamp over the tube. Slide the tube over the port barb and tighten the screw clamp.
7. Attach the amylase dispenser and fill with water to check for leaks.

Note:

Extra tubing and an extra clamp have been provided in case the tubing is damaged during removal. In the event of damage, cut a new length of tubing and also remove the clamp to the “T”. Clamp and install the new tubing and check for leaks.