

To check for leaks at the fittings, do the following:

- 1) Disconnect the solution supply tubes. Leave the Water supply connected.
- 2) Place the bag suspender in place.
- 3) Using hot soapy water, manually fill vessel until the level sensor is covered. (Dish Soap works best).
- 4) Start a Crude Fiber routine. Close the lid.
- 5) Place the open ends of the drain tube and vent line in separate small buckets of water. If you notice "bubbling" in either bucket you have a valve leak.
- 6) After 5-10 minutes, inspect all the areas on the back of the instrument highlighted with a diamond in the image below. Any bubbling indicates a leak and will require a repair.

DELTA

