

Water Heater 120V (A02) Installation

DELTA/DELTAI and A2000/A2000I

Service Procedure # **018**
11-22-22 AES



1. Screw the Vent Tube into the Water Tank Vent Port.
2. Connect either one of the copper lines to the Blue Fitting (Water Supply In) and to your water source.
3. Connect the other tube to the Valve with the Red Fitting (Hot Water Out to the instrument). Turn the water supply on and fill the water heater. Open the Red Fitting to release trapped air. Catch the water in a container. Re-tighten the Red Fitting. Connect the tube to the Instrument's "Water Supply" fitting.
4. Plug the water heater in. If you have a 220v connection, plug your water heater into a Voltage Transformer like the unit shown here (ANKOM part # 6227).

It will take approximately 1/2 hour to reach operating temperature.

5. Refer to the Water Heater owner's manual for additional information.



ANKOM Technology 2052 O'Neil Rd Macedon, NY 14502 PH: (315) 986-8090 FAX: (315) 986-8091