

This procedure explains how to attach the Waste Bottle Assembly (Part# FLEX49) to the ANKOM<sup>FLEX</sup> Analyte Extractor

**1. Screw the cap onto the Waste Reservoir Bottle.**




**2. Connect the shorter opaque tube to the Vessel Vent on the left side of the instrument. Make sure the tube is pushed all the way into the port, then tighten the white nut finger tight.**

Vessel Vent




**3. Connect one of the 10 ft opaque exhaust tubes to the waste assembly exhaust line connection. Make sure the tube is pushed all the way into the exhaust line before hand tightening the nut onto the tube.**

Exhaust line connection

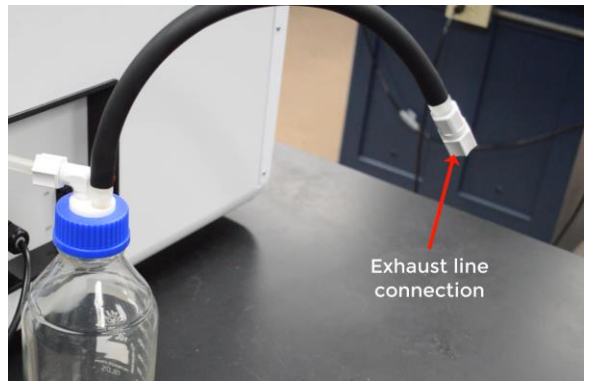


Exhaust tube

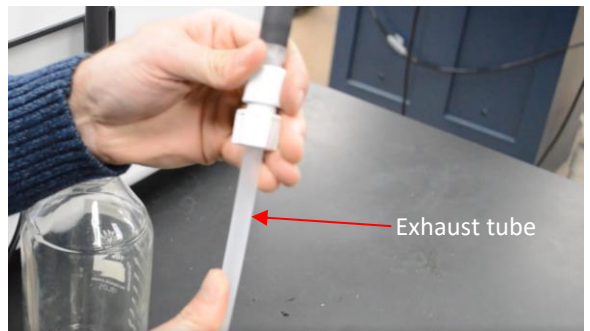

4. Connect the other 10 ft opaque exhaust tube to the globe vent. Slightly unscrew the nut on the fitting to allow the tube to be pushed into the port.

Globe vent

Exhaust tube

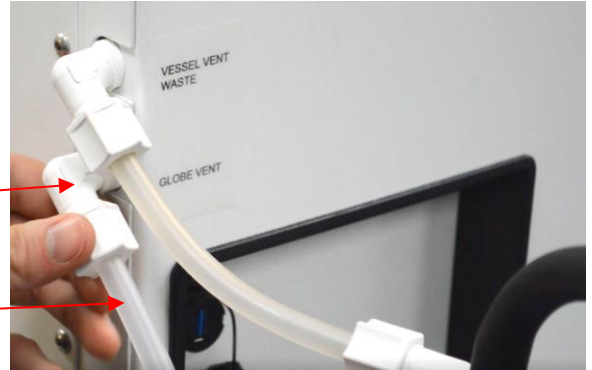

5. Vent both exhaust tubes to exhaust system.

Exhaust Tubes

