


Before assembling the digestion vessels, make sure that each part is clean.

1. **Select the appropriate filter for the analysis. Place the filter (either side up) into the digestion vessel base. Make sure that it is pressed all the way down.**

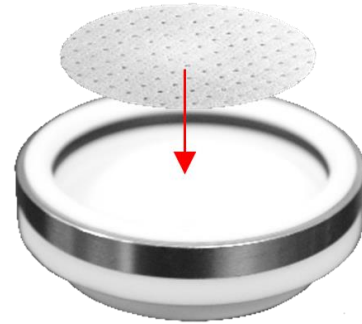

2. **Wet the bottom outside edge of the digestion vessel with water.**



3. **Place the digestion vessel upside-down on the counter. Place the vessel bottom assembly on the digestion vessel and press down firmly until the digestion vessel seats evenly in the vessel bottom assembly.**



IMPORTANT: Make sure the digestion vessel seats evenly and tightly, with the glass all the way down in the digestion vessel base, otherwise the digestion solution will leak.

4. Place the magnetic cross stir-bar into the assembled digestion vessel.



Assembled Digestion Vessel

Digestion Vessel
Glass without Port
(Part # 9362)



Vitamin and Cholesterol
Vessel Filter
(Part # FLEX-VF)
Or
Fat Vessel Filter
(Part # FLEX-FF)



Magnetic Cross Stir-bar
(Part # 9415)



Digestion Vessel Base
(Part # 9407)



Other Digestion Vessel
types:



Digestion Vessel Glass with Port
(Part # FLEX55)