


After the ANKOM<sup>FLEX</sup> method has ended, the round bottom flasks in the recovery oven will contain the isolated analyte.

1. Turn the large black release ring counterclockwise (to the left) to release the round bottom flasks.



2. Remove the round bottom flasks and immediately cover each flask with aluminum foil or a stopper to protect the analyte.

Make sure the covering is secure.

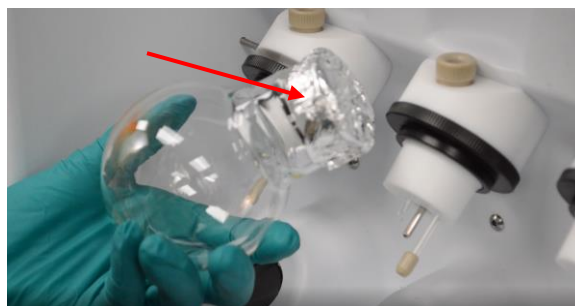

3. Place each covered round bottom flask underneath cold running water to cool the analyte.

Cooling protects the analyte from degradation.



4. Place round bottom flasks in an appropriate stand until ready for post-processing.