

The (4) Vessel Tops in the FLEX Digestion Oven contain O-rings (PN 9231) which provide a seal to the Digestion Vessels and should be replaced if leaking. When a leak occurs, first check that the O-ring is clean and that the mating surface (Vessel) is also clean and undamaged. If the leak persists, then replace the O-ring as follows:

1. Select the 90 Degree Angled Pick PN 9488 from the provided Tool Kit.



2. Gently push the point of the Pick into the rubber O-ring and lift it from the groove. Avoid damaging the groove with the Pick.



3. Clean any residue from the groove with water. Dry the groove and install the new O-ring. Ensure that it is pushed all the way into the groove.

