


Follow the steps below to replace the Exhaust Valve in your ANKOM^{A200/A200I} instrument.




1. Open the back of the instrument.

Lay the A200 on it's front side. Do not put weight on the Controller, Heat or Agitate switches. Use a Phillips screwdriver to remove the four feet and the bottom panel.

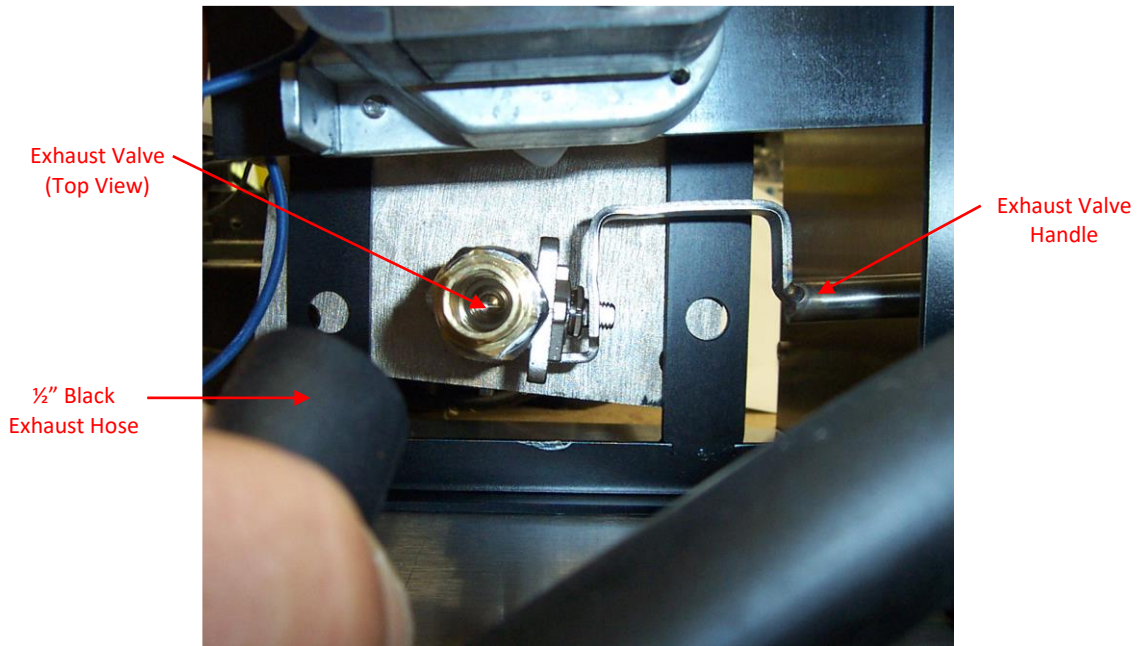

2. Remove the exhaust hose clamp and the 1/2" black exhaust hose.

3. Remove the nut securing the Exhaust Valve Handle to the Exhaust Valve using a 1/2" wrench.

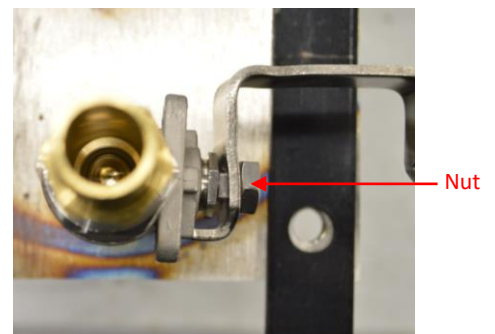

4. Remove the Exhaust Valve by using a large adjustable wrench (turn counterclockwise).

5. Wrap the valve with Teflon Tape.

Use three wraps of Teflon tape on the nipple and install the new valve.



6. Reinstall the Exhaust Valve.

When tightening, orient the opening bolt towards the Exhaust Valve Handle. Tighten the nut holding the Exhaust Valve Handle to the valve.

7. Secure the Exhaust hose with the provided clamp.

8. Re-attach the bottom panel.

With this completed, you are now ready to return your A200/A200I instrument to use.