

The following procedure will instruct you on how to effectively open the F58 bag for sample insertion.

1. **Insert a finger into the slot to open the bag.**




2. **Pinch the sealed sides of the bag to hold it open.**



3. **Press the opposite, pinched sides of the bag together.**




Side View

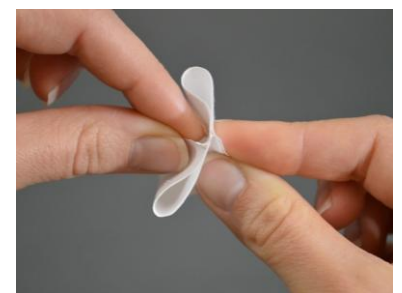

Top View

4. Place the bag on a flat surface with one of the sealed edges face down. Crease the two outside edges to form a V-shape in the material.

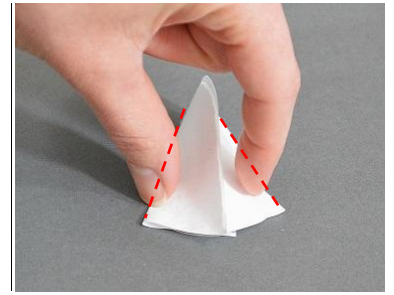

Front View



Side View

5. Place the bag into a bag weigh holder. The improved diamond shape of the bag will help it to remain open more effectively as sample is inserted.

